
TURBULA[®] SYSTEM SCHATZ

Dreidimensionaler Schüttelmischer
Three-dimensional shaker mixer
Mélangeur dynamique tridimensionnel

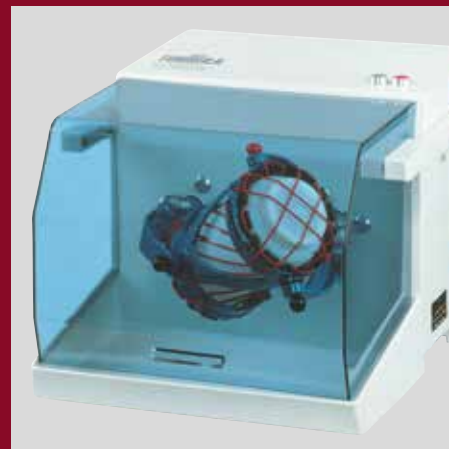Prinzip

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Basic principle

The TURBULA® shaker-mixer is used for the homogeneous mixing of powdery substances with differing specific weights and particle sizes. The product is mixed in its own closed container. It is also possible to mix wet and dry components or different wet components. The production process is hygienic and dust-free, making the TURBULA® easy to clean.

The exceptional efficiency of the TURBULA® shaker-mixer comes from the use of rotation, translation, and inversion according to the Schatz geometric theory.

The mixing container is set into three-dimensional movement that exposes the product to always changing, rhythmically pulsing motion. The results fulfill the highest requirements and are achieved in a minimum of time.

Advantages

- minimum machine maintenance
- dust-free processing
- easy cleaning
- minimum shear force
- no product separation

Application fields

TURBULA® shaker-mixers are used wherever minimum mixing times together with fulfillment of the highest homogeneity standards are required - in all industries, in research and development, and in production plants.

- Diamond tools
- Electrotechnology
- Sintered metals
- Ceramic industry
- Cosmetic industry
- Pharmaceutical industry
- Chemical industry
- Food industry
- Analysis

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TURBULA® T 2 F

The mixing basket can hold any form of container from test tubes up to containers having a maximum volume of 2 liters. The containers are fastened in place by twisted rubber rings. The basket movement is driven by elastic drive belts and an eccentric drive gear. The speed can be varied by adjusting the position of the drive belts on the 5-step pulleys. The speed can also be adjusted by means of an optional frequency converter set up next to the machine.

The TURBULA® T 2 F is used in all industries and application fields, especially in research, development, and analysis.

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TURBULA® T 10 B

The mixing basket can hold standard containers with a maximum volume of 17 liters. Twisted rubber clamping rings allow the use of smaller containers. The movement of the mixing basket is controlled by a silent pendulum chain drive. The speed can be varied by adjusting the position of the round drive belt on the 4-step pulleys.

The TURBULA® T 10 B is often the preferred mixer for pharmaceutical mixing, for mixing sintered metals and ceramic materials, and for many other application fields.

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TURBULA® T 50 A

The mixing basket can hold containers having a maximum volume of 55 liters. Special holders must be mounted for smaller containers. The movement of the mixing basket is controlled by a silent pendulum chain drive. The speed can be varied by adjusting the position of the V-belt on the 5-step pulleys. The mixing basket is brought into loading position by an extra slow speed drive. A hand-cart (optionally delivered with the machine) is used for loading.

The TURBULA® T 50 A is preferred in the pharmaceutical industry, as well as for mixing sintered metals, ceramic materials, and sensitive products.

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A world-wide network with over 40 sales and service points provides support and advice.

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Headquarters of the WAB-GROUP

WAB Machinery (Shenzhen) Co., Ltd.

Unit 102, Floor 1, Plant Building 1
Everbrite Technology Park, Mingqing Road
Longhua District, Shenzhen 518109
China
Tel. +86 755 2813 5108
Fax +86 755 2813 5008
wab-cn@wab-group.com
www.wab-group.com

Willy A. Bachofen GmbH

Liebigstrasse 16
61130 Nidderau-Heldenbergen
Germany
Tel. +49 6187 9394-00
Fax +49 6187 9394-29
wab-de@wab-group.com
www.wab-group.com

WAB India Pvt. Ltd.

Unit No. 1, 2, 3, 4. Plot No. D1/D2
T.T.C. Industrial Area, Turbhe
Navi Mumbai – 400 705
India
Tel. +91 22 613 171-00
Fax +91 22 613 171-10
wab-in@wab-group.com
www.wab-group.com

Willy A. Bachofen S.à.r.l.

3 rue Paul Henri Spaak
68390 Sausheim
France
Tel. +33 389 31 17 00
Fax +33 389 61 89 94
wab-fr@wab-group.com
www.wab-group.com

WAB US Corp.

3 Pearl Court
Suite E
Allendale, NJ 07401
USA
Tel. +1 973 873-9155
wab-us@wab-group.com
www.wab-group.com